
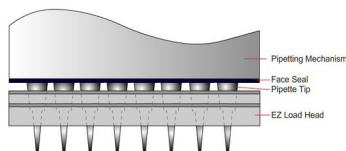


EZ-Load™ Pipettor Heads

The EZ-Load pipettor heads from Apricot Designs simplify the use of disposable tips for multiple channel pipetting by making tip loading and unloading a simple manual process without the requirement for any tools.

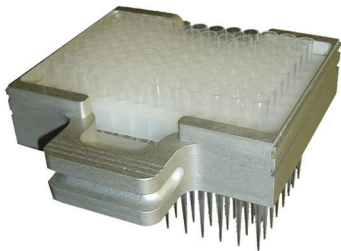

The EZ-Load heads are used in conjunction with specially-made disposable pipette tip racks that allow the transfer of a rack of tips directly from the package into the EZ-Load head. The operator simply drops the rack of tips into the head, places the head in the pipettor, and presses a button to load the head. **Unlike all other tip loading designs, the end of the tip is not forced over a protruding mandrel. No longer is it necessary to apply excess force or “tap” in order to get the tips properly loaded.**

The EZ-Load heads feature a Swing Lock design that allows the tips to be placed into the head without any contact on the bottom part of the tips, thus eliminating cross-contamination due to contact with the head after pipetting. The tops of the pipette tips are sealed against the face seal by the pipettor’s automatic electric loading mechanism. A reliable seal for all tips is provided for optimal pipetting performance.


4 easy steps to load disposable tips and start pipetting!



Remove tips from package rack



Place tips in EZ-Load head

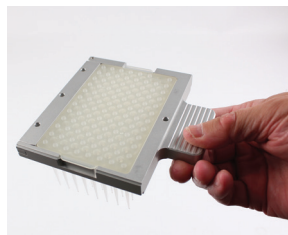

Load head into pipettor



Press the Load/Unload button

Ordering Information

Part Number	Description	Compatible Tips	Compatible Pipettors
096-A-01-EZL-SL-550	EZ-Load Pipetting Head, Swing Lock Design, 96-channel	Apricot 550	PP-96H-M series, i-Pipette96-500 series
096-A-01-EZL-SL-1000	EZ-Load Pipetting Head, Swing Lock Design, 96-channel	Apricot 1000	i-Pipette96-1000 series
096-A-01-EZL-SL	EZ-Load Pipetting Head, Swing Lock Design, 96-channel	Apricot EZL	PP-96L-M series, i-Pipette96-125 series
384-A-01-EZL-SL	EZ-Load Pipetting Head, Swing Lock Design, 384-channel	Apricot EZL	PP-384-M, i-Pipette384 series
384-A-01-EZL-SL-096	EZ-Load Pipetting Head, Swing Lock Design, 96-channel (Offset)	Apricot EZL	PP-384-M, i-Pipette384 series